



Cat. No.	PHC02-20 (Transparent)
Product Description	<p>I.S. CIRCLE WRITER makes a strong circle that will not fall off from the slide while staining procedure and repeated rinsing.</p> <ul style="list-style-type: none">◆ Effective for immunostaining procedures by the Peroxidase-Antiperoxidase(PAP) method, ABC method, B-SA method, immunofluorescence method, ASD method, Enzyme method and Frozen Section method in the histology, cytology, cell culture, biology and microbiology◆ Enable to write any size of circle and pattern on glass slides◆ Proves to be carried out the maximum efficiency with minimum amount of valuable antiserum
Storage	Room Temperature
Usage	<ul style="list-style-type: none">♥ Before use of the pen, hold the pen point up and press it down with fingertip to take out the air in a pen barrel for preventing over flow. It is necessary to perform it by all means.♥ If the glass slide is hydrophile treated, the glass slide should be soaked in a 0.1N hydrochloric acid solution for about 30 seconds and then dry it after washing with water for avoiding fall-off the circle from the slide.♥ Application for <u>Frozen Sections</u>: Before applying the pen, it is necessary to clean up the glass slide by xylene and ethanol or coat with egg-albumin glycerine or poly-L-lysine.